

Cambridge International AS & A Level

BIOLOGY 9700/33

Paper 3 Advanced Practical Skills 1

February/March 2024



This document gives details of how to prepare for and administer the practical exam.


The information in this document and the identity of any materials supplied by Cambridge International are confidential and must NOT reach candidates either directly or indirectly.

The supervisor must complete the report at the end of this document and return it with the scripts.

INSTRUCTIONS

 If you have any queries regarding these confidential instructions, contact Cambridge International stating the centre number, the syllabus and component number and the nature of the query.
 email info@cambridgeinternational.org

phone +44 1223 553554

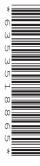

General information about practical exams

Centres must follow the guidance on science practical exams given in the Cambridge Handbook.

Safety

Supervisors must follow national and local regulations relating to safety and first aid.

Only those procedures described in the question paper should be attempted.

Supervisors must inform candidates that materials and apparatus used in the exam should be treated with caution. Suitable eye protection should be used where necessary.

The following hazard codes are used in these confidential instructions, where relevant:

CcorrosiveMHmoderate hazardHHhealth hazardTacutely toxicFflammableOoxidising

N hazardous to the aquatic environment

Hazard data sheets relating to substances used in this exam should be available from your chemical supplier.

Before the exam

- The packets containing the question papers must **not** be opened before the exam.
- It is assumed that standard school laboratory facilities, as indicated in the *Guide to Planning Practical Science*, will be available.
- Spare materials and apparatus for the tasks set must be available for candidates, if required.

During the exam

- It must be made clear to candidates at the start of the exam that they may request spare materials and apparatus for the tasks set.
- Where specified, the supervisor must perform the experiments and record the results as instructed.
 This must be done out of sight of the candidates, using the same materials and apparatus as the candidates.
- Any assistance provided to candidates must be recorded in the supervisor's report.
- If any materials or apparatus need to be replaced, for example, in the event of breakage or loss, this must be recorded in the supervisor's report.

After the exam

- The supervisor must complete a report for each practical session held and each laboratory used.
- Each packet of scripts returned to Cambridge International must contain the following items:
 - the scripts of the candidates specified on the bar code label provided
 - the supervisor's results relevant to these candidates
 - the supervisor's reports relevant to these candidates
 - seating plans for each practical session, referring to each candidate by candidate number
 - the attendance register.

Specific information for this practical exam

During the exam, the supervisor or other competent biologist (**not** the invigilator) should obtain the results needed for the supervisor's report by following the relevant steps in the question paper. The results should be recorded in the supervisor's report.

Organisation of the exam

- All candidates must have access to the materials required for Question 1 throughout the whole period of the exam.
- Half of the candidates will have access to the microscope and slide for a maximum time of one hour from the start of the exam. These candidates should start with Question 2. After one hour, or sooner if candidates have finished Question 2, they should move on to Question 1.
- For Question 2, two candidates are **not** permitted to share the same microscope and slide at the same time.
- The other candidates should start with Question 1. After one hour, these candidates should be given access to the microscope and slide. They should then move on to Question 2 as soon as they are ready.
- Candidates will only have access to the microscope and slide for one hour. They should be
 advised that they can answer any part of the exam paper not requiring the microscope and slide
 throughout the whole period of the exam.
- Access arrangements to microscopes and slides, including instructions on which question to start with and timings, must be explained to candidates before the start of the exam.

Materials to be supplied by Cambridge International

Slide P1

On receipt of the slides, check that they are labelled **P1** and that no slides are broken. The slides should **not** be viewed in advance of the exam. The material on the slides is confidential and must **not** be disclosed to candidates.

The number of slides supplied by Cambridge International will be equal to half the candidate entry.

Return of slides to Cambridge International

Immediately after the exam, the slides must be:

• returned to Cambridge International in the boxes in which they were received, using the self-adhesive label supplied. The slides must **not** be included in the packet of scripts.

or

purchased using the order form enclosed with the slides, which should be completed and returned
to Cambridge International. The order form must **not** be included in the packet of scripts. Slides
and boxes will be charged at the rate of £3.25 per slide plus £1 per box.

If the slides are not returned or purchased by the deadline stated on the order form, the charge will be £3.75 per slide plus £1 per box.

Materials and apparatus for Question 1

Each candidate will need:

0.5% glucose solution in a beaker, labelled G1, provided at room temperature (see Preparation of materials) 0.25% glucose solution in a beaker, labelled H1, provided at room temperature (see Preparation of materials) 10.25% glucose solution in a beaker, labelled H1, provided at room temperature (see Preparation of materials) 10.25% glucose solution in a beaker, labelled H1, provided at room temperature 10.25% glucose solution of materials) 10.25% glucose solution in a beaker, labelled H1, provided at room temperature 21.25 cm³ 22.26 cm³ 23.26 cm³ 24.27 cm³ 25.27 cm³ 26.28 eakers, capacity 50–100 cm³ 27.28 eakers, capacity 50–100 cm³ 28.29 eakers, capacity 50–100 cm³ 29.30 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath 10.25 cm² 10.25 cm³ 10	materials and apparatus for each candidate	quantity	1
(see Preparation of materials) 10 cm³ 0.25% glucose solution in a beaker, labelled H1, provided at room temperature (see Preparation of materials) at least 10 cm³ Distilled water in a beaker, labelled W, provided at room temperature at least 150 cm³ [MH][N] Benedict's solution in a beaker, labelled Benedict's, provided at room temperature at least 25 cm³ 10 cm³ syringes 2 5 cm³ syringes 2 8 eakers, capacity 50–100 cm³ 5 Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath 1 Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels 8 Glass marker pen (permanent) 1 Glass rod 1 Stop-clock or timer showing seconds 1	1.0% glucose solution in a beaker, labelled R , provided at room temperature (see Preparation of materials)		
(see Preparation of materials) Distilled water in a beaker, labelled W, provided at room temperature [MH][N] Benedict's solution in a beaker, labelled Benedict's, provided at room at least temperature [MH][N] Benedict's solution in a beaker, labelled Benedict's, provided at room at least temperature 10 cm³ syringes 2 Seakers, capacity 50–100 cm³ Seakers, capacity 50–100 cm³ Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Stop-clock or timer showing seconds	0.5% glucose solution in a beaker, labelled G1 , provided at room temperature (see Preparation of materials)		
Distilled water in a beaker, labelled W, provided at room temperature 150 cm ³ MH][N] Benedict's solution in a beaker, labelled Benedict's, provided at room temperature 25 cm ³ 10 cm ³ syringes 2 5 cm ³ syringes 2 Beakers, capacity 50–100 cm ³ 5 Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm ³ , suitable for heating as a water-bath, with approximately 200 cm ³ of tap water at 40–50 °C, labelled water-bath 7 Test-tubes, small, capacity 20–30 cm ³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm ³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm ³ , labelled For waste 1 Paper towels 8 Glass marker pen (permanent) 1 Glass rod 1 Stop-clock or timer showing seconds 1	0.25% glucose solution in a beaker, labelled H1 , provided at room temperature (see Preparation of materials)		
temperature 25 cm³ 10 cm³ syringes 2 5 cm³ syringes 2 Beakers, capacity 50–100 cm³ 5 Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath 7 Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more 1 Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels 8 Glass marker pen (permanent) 1 Stop-clock or timer showing seconds 1	Distilled water in a beaker, labelled W , provided at room temperature		
5 cm³ syringes Beakers, capacity 50–100 cm³ 5 Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Stop-clock or timer showing seconds	[MH][N] Benedict's solution in a beaker, labelled Benedict's , provided at room temperature		
Beakers, capacity 50–100 cm³ Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Stop-clock or timer showing seconds	10 cm ³ syringes	2	
Bunsen burner, tripod, gauze, bench mat and lighter or matches 1 Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ 7 Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Stop-clock or timer showing seconds	5 cm ³ syringes	2	
Beaker, capacity approximately 400 cm³, suitable for heating as a water-bath, with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Stop-clock or timer showing seconds	Beakers, capacity 50–100 cm ³	5	
with approximately 200 cm³ of tap water at 40–50 °C, labelled water-bath Test-tubes, small, capacity 20–30 cm³ Test-tube rack(s) to hold 7 small test-tubes 1 or more Test-tube holder to hold hot test-tubes 1 Container, with approximately 200 cm³ of tap water, labelled For washing 1 Container, capacity approximately 400 cm³, labelled For waste 1 Paper towels Glass marker pen (permanent) 1 Glass rod 1 Stop-clock or timer showing seconds	Bunsen burner, tripod, gauze, bench mat and lighter or matches	1	
Test-tube rack(s) to hold 7 small test-tubes Test-tube holder to hold hot test-tubes Container, with approximately 200 cm³ of tap water, labelled For washing Container, capacity approximately 400 cm³, labelled For waste Paper towels Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds 1 or more 1 or more	Beaker, capacity approximately 400 cm ³ , suitable for heating as a water-bath, with approximately 200 cm ³ of tap water at 40–50 °C, labelled water-bath	1	
Test-tube holder to hold hot test-tubes Container, with approximately 200 cm³ of tap water, labelled For washing Container, capacity approximately 400 cm³, labelled For waste Paper towels Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds	Test-tubes, small, capacity 20–30 cm ³	7	
Container, with approximately 200 cm³ of tap water, labelled For washing Container, capacity approximately 400 cm³, labelled For waste Paper towels Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds 1	Test-tube rack(s) to hold 7 small test-tubes	1 or more	
Container, capacity approximately 400 cm ³ , labelled For waste Paper towels Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds 1 1 1 1 1	Test-tube holder to hold hot test-tubes	1	
Paper towels Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds 8 1	Container, with approximately 200 cm ³ of tap water, labelled For washing	1	
Glass marker pen (permanent) Glass rod Stop-clock or timer showing seconds 1	Container, capacity approximately 400 cm ³ , labelled For waste	1	
Glass rod 1 Stop-clock or timer showing seconds 1	Paper towels	8	
Stop-clock or timer showing seconds 1	Glass marker pen (permanent)	1	
3 2 2 2 3 3 3 3 3 3 3 3 3 3 3 3 3 3 3 3	Glass rod	1	
Suitable eye protection 1	Stop-clock or timer showing seconds	1	
	Suitable eye protection	1	

Preparation of materials

All the materials may be prepared the day before the exam.

Solutions R, G1 and H1 should be kept overnight in a covered container in a refrigerator.

All solutions and the distilled water, **W**, should be at room temperature before the start of the exam.

• R, 1.0% glucose solution

This is prepared by putting 1g of glucose into 50 cm³ of distilled water. Make up to 100 cm³ with distilled water. Mix well.

• **G1**, 0.5% glucose solution

This is prepared by putting 0.5 g of glucose into 50 cm³ of distilled water. Make up to 100 cm³ with distilled water. Mix well.

• **H1**, 0.25% glucose solution

This is prepared by putting 0.25 g of glucose into 50 cm³ of distilled water. Make up to 100 cm³ with distilled water. Mix well.

Materials and apparatus for Question 2

Each candidate will need:

materials and apparatus for each candidate	quantity	✓
Microscope with: an eyepiece lens, ×10 magnification a low-power objective lens, ×10 magnification a high-power objective lens, ×40 magnification 	1 between 2	
Slide P1	1 between 2	
Ruler in mm	1	

Preparation of materials

Microscope

Any lenses which are **not** ×10 or ×40 should be removed or replaced.

For each candidate:

- the microscope must be set up on low power
- the slide must **not** be on the stage of the microscope.

Permission to reproduce items where third-party owned material protected by copyright is included has been sought and cleared where possible. Every reasonable effort has been made by the publisher (UCLES) to trace copyright holders, but if any items requiring clearance have unwittingly been included, the publisher will be pleased to make amends at the earliest possible opportunity.

To avoid the issue of disclosure of answer-related information to candidates, all copyright acknowledgements are reproduced online in the Cambridge Assessment International Education Copyright Acknowledgements Booklet. This is produced for each series of examinations and is freely available to download at www.cambridgeinternational.org after the live examination series.

Cambridge Assessment International Education is part of Cambridge Assessment. Cambridge Assessment is the brand name of the University of Cambridge Local Examinations Syndicate (UCLES), which is a department of the University of Cambridge.

© UCLES 2024

3

Supervisor's report

Syllabus and component number	9	7	0	0	/	3

Centre number

Centre name

Time of the practical session

Laboratory name/number

Give details of any difficulties experienced by the centre or by candidates (include the relevant candidate names and candidate numbers).

You must include:

- any difficulties experienced by the centre in the preparation of materials
- any difficulties experienced by candidates, e.g. due to faulty materials or apparatus
- any specific assistance given to candidates.

Temperature of exam room	°C
Results for Question 1(a)(ii)	

Results for Question 1(a)(iv)

Declaration

- 1 Each packet that I am returning to Cambridge International contains all of the following items:
 - the scripts of the candidates specified on the bar code label provided
 - the supervisor's results relevant to these candidates
 - the supervisor's reports relevant to these candidates
 - seating plans for each practical session, referring to each candidate by candidate number
 - the attendance register.
- 2 Where the practical exam has taken place in more than one practical session, I have clearly labelled the supervisor's results, supervisor's reports and seating plans with the time and laboratory name/number for each practical session.
- 3 I have included details of difficulties relating to each practical session experienced by the centre or by candidates.
- 4 I have reported any other adverse circumstances affecting candidates, e.g. illness, bereavement or temporary injury, directly to Cambridge International on a *special consideration form*.

Signed	. (supervisor)
Name (in block capitals)	

© UCLES 2024 9700/33/CI/F/M/24